

# MouseOx - Collar clip Sensor for awake subjects

Collar clip sensor to collect vital sign measurements, with MouseOx Plus, on non-anesthetized subjects

**Product codes:**

Reference: 01502

**Product description:**

These sensors are in the form of a clip which is applied on the neck of the animal.

They are available in 6 different sizes from 1 to over 500 g. Each sensor is delivered with its blank equivalent, without sensor, so the animal can get used to wearing it before starting the measurements.

**Collar size** Unshaven mouse Shaven mouse Rat **XS** 1-15 g 1-20 g N/A **S** > 15 g > 20 g 5 - 30 g **M** N/A N/A 31 - 100 g **L** N/A N/A 101 - 300 g **XL** N/A N/A 301 - 500 g **2XL** N/A N/A > 500 g

These sensors are designed to provide the **highest quality measurements**. To assure the highest level of accuracy it is recommended that you replace each sensor after **100 hours of use**. Each sensor is shipped with a "Handle Support" in place across the finger tabs to prevent plastic deformation that can accumulate over long periods of time. When not in use always store the sensors with the handle clamp in place.

The CollarClip, which is colored blue, is to be applied **from the back of the neck** of an animal, and can be used in both awake and anesthetized applications.

For awake subject, you need to have the **MouseOx Plus-Premium software**.

### Product features:

### Available models

Description: XS CollarClip (carotid artery) - Conscious Unshaved Mouse 1-15 g, Shaved Mouse 1-20 g (015020), XS CollarClip (carotid artery) - Conscious Unshaved Mouse 1-15 g, Shaved Mouse 1-20 g (015020-5), S CollarClip (carotid artery) - Conscious Unshaved Mouse + 15 g, All Mice + 20 g, Rat 5-30 g (015021), S CollarClip (carotid artery) - Conscious Unshaved Mouse + 15 g, All Mice + 20 g, Rat 5-30 g (015021-5), M CollarClip (carotid artery) - Conscious Rat 31-100 g (015022), M CollarClip (carotid artery) - Conscious Rat 31-100 g (015022-5), L CollarClip (carotid artery) - Conscious Rat 101-300 g (015023), L CollarClip (carotid artery) - Conscious Rat 101-300 g (015023-5), XL CollarClip (carotid artery) - Conscious Rat 301-500 g (015024), XL CollarClip (carotid artery) - Conscious Rat 301-500 g (015024-5), 2XL CollarClip (carotid artery) - Conscious Rat + 500 g (015025), 2XL CollarClip (carotid artery) - Conscious Rat + 500 g (015025-5)

Packing: single unit (015020), pack of 5 pce (015020-5), single unit (015021), pack of 5 pce (015021-5), single unit (015022), pack of 5 pce (015022-5), single unit (015023), pack of 5 pce (015023-5), single unit (015024), pack of 5 pce (015024-5), single unit (015025), pack of 5 pce (015025-5)

# LAB